Certifications by ---

- CE Marked Products.
- ISO 9001 Certified.
- PL Insurance.

◇ English Page : P. 1 ~ 18

◇ Korean Page : P. 19 ~ 36

User Manual

(Version 3.1.2)

written for

WUC-A01 / -A02 / -A03 / -A06 / -A10 / -A22

" WUC-A10 " Analog Ultrasonic Cleaners

Sales & Basic A/S :
By the Nearest Official Distributors.

EUROPEAN DECLARATION OF CONFORMITY

We, DAIHAN Scientific Co., Ltd.,
326, Sinpyeongseokhwa-ro, jijeong-myeon, Wonju-si, Gangwon-do, Korea
Tel: 82-2-967-5235, Fax: 82-2-967-5258
declare under our sole responsibility that the product

Ultrasonic Cleaner **WUC-A01, WUC-A02, WUC-A03,** **WUC-A06, WUC-A10, WUC-A22**

to which this declaration relates is in conformity
with the following standard or other normative document

EN 61326, EN 61000-3-2, EN 61000-3-3
EN 61010-1, EN 61010-2

following the provisions of the directives:

Low-Voltage Directive 2014/35/EU **EMC Directive 2014/30/EU**

EC Directive 2006/42/EC
RoHS - Directive 2011/65 EU

Signed: _____ (Shu, Hyun Jeong)

Date: 01, Januaary 2023

Authority: President / CEO

For

DAIHAN Scientific Co., Ltd.
326, Sinpyeongseokhwa-ro, jijeong-myeon, Wonju-si, Gangwon-do, Korea
Tel: 82-2-967-5235, Fax: 82-2-967-5258

Table of Contents

	Page
1. General	
1.1 Precautions Before Use	4
1.2 Product Introduction	5
1.3 Product Configuration	5
2. Installation	
2.1 Product Installation	6
2.2 Product Specifications	7
3. Product Usage	
3.1 Product Usage	8
3.2 Names of Each Part	11
3.3 Usage Instructions	13
4. Products Management Methods	
4.1 Accessories	16
4.2 Maintenance and Management	16
4.3 Troubleshooting	17
5. Product Assurance	
5.1 Quality Assurance	18

1. General

1

1.1 Precautions Before Use

Thank you for purchasing the Analog Ultrasonic Cleaners.

This manual describes the performance and usage instructions of the product, and the precautions in the handling. Please carefully read this manual, before using the product. The following warnings should be strictly adhered to: (The following warning sign is marked in the part that is required user attention, so please safely use the product after being well-acquainted with the meaning.)

	Warning Failure to follow the instructions may result in injury.
	Caution Failure to follow the instructions may result in product failure.
	Notification Do not disassemble or modify the product or use it for purpose other than its intended function.
	Caution: High Temperature High-temperature heating, pose a risk of burns.
	Danger of Explosion If volatile, explosive chemicals are used, there is a danger of explosion.
	Proximity Caution Keep clear of the product during use.
	This symbol indicates Protective earth (ground).
	This symbol indicates power "OFF" on the power on/off switch.
	This symbol indicates power "ON" on the power on/off switch.
	This symbol indicates the connection point of the neutral conductor on the installed unit.
	This symbol indicates alternating current Voltage.

1. General

1.2 Product introduction

Ultrasonic Cleaner uses for cleaning with the various kinds of dirt like finger marks or oils on the surface located in the irregular narrow spaces at the corner. Built in Heat-on mode to maximize the cleaning efficiency. This product was developed and manufactured considering the best possible ease for use and safety, and has the following features.

1. Can expect the cleansing effect due to heating equipped with a temperature control.
2. 40KHz ultrasonic frequency for Excellent washing effect supports.
3. Stainless steel Bath and Lid is to minimize corrosion.
4. offers a wide range of experimental environment equipped with a timer.
5. CE Certified Products.
6. PL(Product Liability) Insurance

1. General

1.3 Product Configuration

Component Photo	Component Name	No. of Configuration
	Analog Ultrasonic Cleaners " WUC-A01 "	1 Unit
	Analog Ultrasonic Cleaners " WUC-A02 "	1 Unit
	Analog Ultrasonic Cleaners " WUC-A03 / -A06 "	1 Unit
	Analog Ultrasonic Cleaners " WUC-A10 / -A22 "	1 Unit
	Flat Lid	1 ea
	Spare Fuse	2 ea

-Including a copy of the Operation Manual

2. Installation

2

2.1 Product Installation

1. Please do not lay the product on the floor or subject it to shock. It can cause damage inside the product.
2. Please install the product in a hard, flat, and stable place.
3. Please do not install the product in a place where it is exposed to direct sunlight, or in the dangerous places.
4. To minimize the risk of short circuits, do not install the product (in the) places where it will be subjected to the inflow of moisture or organic solvents on the inside of the body.
5. This product was manufactured for normal operation at a rated voltage, so please check the voltage status before installing the product.
6. When installing and using the product, please certainly use the power cord provided with the products.
7. Please be careful so that moisture, organic solvent, dust, corrosive and gas, etc. do not enter into this product.
8. Please do not expose this product in the place where there is direct heat sources.

* Caution

- The warnings sticker is affixed to the product. Please read.

Example) WUC-A10

* WUC-A01

- Do not make water level below 3 c m
- Do not operate continuously more than 5 minutes
- Keep 5 minutes break time when after 5 minutes run continuously.

* WUC-A02H/A03

- Do not make water level below 6 c m
- Do not operate continuously more than 30 minutes
- Keep 10 minutes break time when after 30 minutes run continuously.

* WUC-A06/A10/A22

- Do not make water level below 10 cm
- Do not operate continuously more than 40 minutes
- Keep 10 minutes break time when after 40 minutes run continuously.

* Common Point

- Turn off the heater power if the unit is not going to be used.
- When make water level below standard, it may cause damage of heater.
- Do not use acid detergent solution(Chlorine line).

2. Installation

2.2 Product Specifications

Models	WUC-A01	WUC-A02	WUC-A03	WUC-A06	WUC-A10	WUC-A22
Capacity	1.2 Lit	1.8 Lit	3.3 Lit	6 Lit	10 Lit	22 Lit
Frequency	40 KHz					
HF-Peak Out	72 W	75 W	124 W	168 W	265 W	379 W
Heating Temp.	65°C Fixed		45°C to 85°C			
Heating Power	86 W	86 W	172 W	172 W	258 W	275 W
Timer	1~5 min	5~30 min	0~40 min			
Drain Valve	None				1/2" NPT	
Tank Size inner (w x d x h) (mm)	150X135X60	150X135X100	240X140X100	290X150X150	290X240X150	500X300X150
Weight	1.7 kg	2.7 kg	4.2 kg	5.7 kg	7.8 kg	12 kg
Packing Dimension (w x d x h) (mm)	235X205X215	228X205X240	310X225X298	380X240X360	450X335X380	650X450X400
Gross Weight	2.1 kg	3.1 kg	5 kg	6.7 kg	9.4 kg	14.3 kg
Power Consumption	115 W	115 W	230 W	280 W	528 W	739 W
Power Supply	1 Phase, AC 120V, 60Hz or AC 230V, 50/60Hz					

* Other Specifications are available upon Customer's Request.

Product Liability (PL) Insurance Admission

DAIHAN Scientific (Co., Ltd.) is a member of Product Liability Insurance (PL Insurance), so that its customers can trust our products, (believe and use its products,) by preparing against the accident which can be occurred during using the products by self-production.

[Contents]

Insurance Type: Product Liability Insurance

Publishing Office: Machinery Insurance Cooperative

Limit of Indemnity: 100 million won for Persons and Objects

Secured Area: Worldwide (excluding North America)

Emergency Contact: **DAIHAN Scientific (Co., Ltd.) (TEL: 080-008-3000, FAX: 033-737-7575)**
(TEL: 02-369-8516, FAX: 02-369-8521)

3. Product Usage

3

3.1 Precautions for Use

	<p>Please be familiar with the user manual before using this product, and operation should be performed only by qualified personnel</p>
	<p>This product should be connected to a appropriate power source that it is generally used. Please use a lower power source than the regulation load permitted (AC 120V~, 60Hz or AC 230V~, 50/60Hz, Single-phase).</p>
	<p>Place the product on a level, horizontal surface to support the weight of the product without vibration. Install the product in a place where the floor is hard and flat. Failure to do so can cause accidents. by negligence.</p>

● **Warnings**

- **Please observe the following precautions.**

	<p>This product should be operated in a room with temperature maintain between 5℃ and 40℃. Long-term use in an environment that exceeds the allowable range may cause a malfunction.</p>
	<p>Please do not damage or distress the power cord. (Please do not hurt on the power cord, or process it, or force bend it, or pull it, and tie it.) Also, if heavy objects are placed over the power cord, or the power cord is inserted in between objects, the power cord can be damaged and may cause fire or electric shock.</p>
	<p>Please do not directly spray water on the product or wash it with water, to avoid short circuits or electric shock.</p>
	<p>Please do not touch or manipulate the electrical parts such as the power plug with wet hands electric shock.</p>
	<p>Please do not operate the product by manipulation of the plug to avoid electric shock or short circuits</p>
	<p>Do not install the product where it can get damp or in a dusty environment. It may cause overheating or short circuit. – Pollution Degree 2</p>
	<p>Do not install the product in a location at an altitude of over 2,000 meters (6,562 feet).</p>
	<p>Please do not operate in the condition that explosive or flammable chemicals (alcohol, benzene and other flammable solvent) are put inside the Bath.</p>

	<p>If you raise your hand when operating at high temperatures, you may get burned. Wear safety equipment before using it.</p>
	<p>The temperature of the regulator SAFETY should be set up 20°C higher than the operating temperature. If the regulator SAFETY is set to 0, the heater does not operate.</p> <p>If unexpected sound, smell or smog is generated, pull the main plug out from the wall socket and contact the technical assets of DAIHAN Scientific</p> <p>If irradiated or contaminated samples are placed in the chamber, the warranty is voided. Do not use this product to sterilize or disinfect objects or samples.</p>
	<p>Do not use outdoors avoid overheating or electrical faults. The use in the place where the sun continuously shines and can get rained on it is prohibited. It may cause overheating or electrical fault of the product.</p>
	<p>Please do not use flammable or volatile products in the vicinity of the product to minimize the risk of fire. Please do not use the flammable spray such as lacquers and paint, or volatile, flammable chemicals and its similar product, in the proximity of the product. They may cause the ignition due to the spark of switch.</p>
	<p>Please avoid the disassembly or modification of this product by a non-authorized person to avoid serious electrical problems and to retain the warranty service of DAIHAN Scientific Co., Ltd.</p>
	<p>This product must be connected to a properly grounded power supply.</p>
	<p>Please regularly check whether dust or dirt are on the power plug and completely connect the plug. If the plug is covered with dust, or the connection is not complete, it will cause an electrical shock or fire.</p> <p>Please separate the power plug from the outlet for safety when the product is unused for long periods of time to avoid overheating or ignition due to dust.</p> <p>When disconnecting the power plug, it must be grasped by the hands and then pulled. Grasping only the cord and then disconnecting it may cause overheating or ignition because of internal lines being disconnected.</p> <p>Insert the plug securely into an appropriate wall socket. Loose connections result in excessive heat generation to the device and may cause arcing to occur at the connection.</p> <p>If unusual noise or signs of abnormal behavior are found during use of this product, please immediately contact the supplier or DAIHAN Scientific (Co., Ltd.) after turning off the power supply and disconnecting the power line.</p> <p>When moving the product, please move it so that the power cord and the product</p>

	are not damaged after disconnecting the power plug from the electrical outlet, in order to avoid the electrical shock or the ignition due to the damage to the cord.
	Please maintain the clean condition of this product by removing the foreign matter while gently wiping with the neutral detergent and a dry cloth after checking whether the foreign matter present or not, before and after using this product.
	The fluctuation of main power voltage should be lower than +/-10% of the nominal voltage..
	The maximum relative humidity is 80% at 31°C and 50% at 40°C .
	If a product was transported or stored in high humidity condition, please check packaging condition before use. If the carton or wooden box is waterlogged, please contact the deliverer or our technical service engineer. Do not operate the unit before checking with an engineer, otherwise it may cause an electrical shock or a fire.
	The Sound Level of the buzzer is maximum 60dB in 1 meter distance when an error is detected.
	The device should be set at a well ventilated place. Be sure the holes on the side or rear surface of the body are blocked by a wall or an object. And do not operate the device at a poor ventilated place or use closed. It may cause a fire or damage of the device as not to release heat or insufficient ventilation may result.
	If you don't use the device for a long time, please drain all of water from bath and then clean up.
	Never insert the explosive or flammable chemicals inside the Bath. There is a fire danger.
	Please apply only distilled water to the Bath. If you don't, it may cause roast by ferrum which was in tap water.

※ All of the circuits, electrical or mechanical parts and configurations used in this product are the technical assets of DAIHAN Scientific (Co., Ltd.). Only technicians of the technical support team at DAIHAN Scientific (Co., Ltd.) or technicians approved through the education from DAIHAN Scientific should be allowed to repair this product.

3. Product Usage

3.2 Names of Each Part


① WUC-A01

NO	Name	Description
①	Heater On Switch	Running or stopping the heating motion
②	Ultrasonic On & Timer Switch	Run ultrasonic operation and set timer time.
③	Power On Switch	Turn it on/off.

< Figure 3.2.1> The appearance of WUC-A01

② WUC-A02

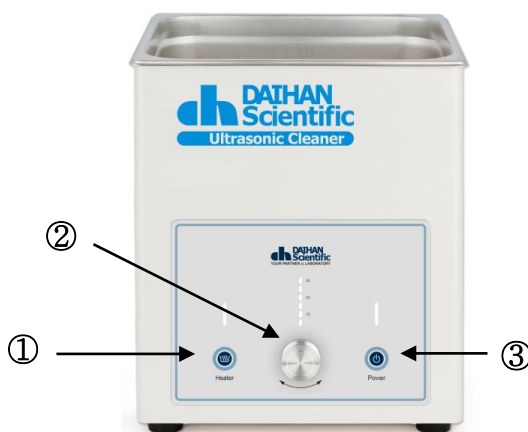
NO	Name	Description
①	Heater On Switch	Running or stopping the heating motion
②	Ultrasonic On & Timer Knob	Run ultrasonic operation and set timer time.
③	Power On Switch	Turn it on/off.

< Figure 3.2.2> The appearance of WUC-A02

③ WUC-A03, A06

④ WUC-A10, A22

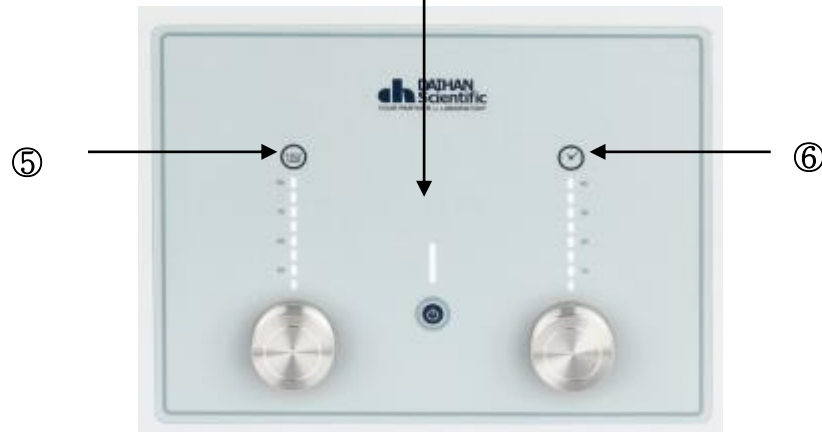NO	Name	Description
①	Heater On & Temp. Knob	Running the heating operation and setting the temperature setting.
②	Power On Switch	Run ultrasonic operation and set timer time.
③	Ultrasonic On & Timer Knob	Turn it on/off.
④	Drain Valve	Drain the solution in the water tank.
⑤	Heter On & Temp. Lamp	Display during the heating motion.
⑥	Ultrasonic On & Timer Lamp	Display during ultrasound
⑦	Power On Lamp	Turn it on/off.

< Figure 3.2.3> The appearance of WUC-A03, A06, A10, A22

3. Product Usage

3.3 Usage Instructions

* Caution

- Ultrasonic Cleaner is 2 types are available.
 - 1) Direct type : Clean the targets in the tube directly.
 - 2) Indirect type : Clean the targets in the tube after clean outside beaker or cup.
(This type are used for acid or corrosion materials.)
- Do not operate over 30min.
- Do not operate without liquid in the Bath.
- Please use water or water and washing solution mix of accordance with the washing object.

* Usage Instructions

- Please put in the bath to water or water and washing solution mix of accordance with the washing object. (About 2/3)
- (Please search for "concentrated cleaner" ALL for LAB.com or www.DAIHAN-Sci.com search a variety of items than the cleaners.)
- Put the stuff to be clean in a bath.
- Connect the power cord to the proper power socket and Turn on the Main Power Switch.
- The operating of product according to the method used by the model below.
- After the completion of cleaning, take out the cleaned materials. Rinse them carefully to remove all the remained cleaning liquid on them, and wipe out with towel.

① WUC-A01

< Figure 3.3.1> Usage Instructions of WUC-A01

- When performing ultrasonic cleaning, turn on the power of the ultrasonic cleaner by pressing the ultrasonic operation switch located in the front control unit of the product for 3 seconds. After cleaning is completed, press the ultrasonic operation switch to turn it off. The

power will be automatically turned off after 3 hours due to the safety function.

- When performing the thermal operation, the heater operates by pressing the thermal operation switch located in the front control part of the product. After cleaning is completed, press the warm operation switch to turn off the heater operation.
- Keep the temperature of the cleaning liquid 65 °C Heater operation.
- It operates when the ultrasound operation switch is pressed, and it operates for 1 minute when pressed once, 3 minutes when pressed twice, and 5 minutes when pressed three times, and the lamp turns on. Even if the power is off, you will remember the last move you used.

② WUC-A02

< Figure 3.3.2> Usage Instructions of WUC-A02

- Press the ultrasonic operation switch for 3 seconds to power on the ultrasonic cleaner. The power will be automatically turned off after 3 hours due to the safety function..
- It works when you press the ultrasound operation Knob, and when you turn the Knob, the timer increases or decreases. You can set it up to 30 minutes every 5 minutes.
- When performing the thermal operation, the heater operates by pressing the thermal operation switch located in the front control part of the product. After cleaning is completed, the heater operation stops when the warm operation switch is turned off..
- Keep the temperature of the cleaning liquid 65 °C Heater operation.

③ WUC-A03H, A06H, A10H, A22

< Figure 3.3.3> Usage Instructions of WUC-A03, A06, A10, A22

- Press the power switch for 3 seconds to power on the ultrasonic cleaner. The power will be automatically turned off after 3 hours due to the safety function.
- It works when you press the ultrasound operation Knob, and when you turn the Knob, the timer increases or decreases. You can set it up to 40 minutes every 5 minutes.
- The heater operates when the thermostat is pressed during the warm-up operation. If you turn the knob, you can adjust it from 45°C to 80°C, and when stopping the operation, press the regulator to turn it off.

***notification**

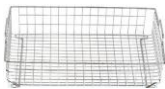
- **This product can set temperature and heating, but does not guarantee temperature accuracy. This is because the sonicator heats in the product.**
- **The heating function is a function for improving cleaning power and does not cover temperature accuracy.**

4. Products Management

4

4.1 Accessories

*** Ordering Information**

Image	Ordering Number	Information
	DH. WUC.B01	Basket wire, "B01", for 1.2Lit, 144 x 131 x h54mm
	DH. WUC.B02	Basket, Wire, "B02", for 1.8Lit, 144 X 131 X h92mm
	DH.WUC.B03	Basket, Wire, "B03", for 3.3Lit, 224 X 128 X h83mm
	DH.WUC.B06	Basket, Wire, "B06", for 6Lit, 267 X 136 X h126mm
	DH.WUC.B10	Basket, Wire, "B10", for 10Lit, 276 X 218 X h132mm
	DH.WUC.B22	Basket, Wire, "B22", for 22Lit, 467 X 275 X h131mm
	DH.WUC.B100	Fine Mesh Strainer, "B100", for Ultrasonic Cleaner, ø100mm

4. Products Management Method

4.2 Maintenance and Management

Analog Ultrasonic cleaners were created in consideration of various situations that may occur during operation. However, in order to use it for a long time with stable performance, please follow the following.

1. Be sure to use rated power for AC 120 or AC 230V, 50/60Hz, and Single phase.
2. Operate indoors without moisture as much as possible, and maintain the indoor temperature at 5°C to 40°C.
3. If not used for a long time, remove the power plug, dry it, and wrap it and store it.
4. If parts need to be replaced due to a failure during use, only genuine products provided by our company should be used.
5. Failure due to use beyond normal limits cannot be maintained.
6. Please use only the prescribed power.
7. Do not touch the power plug and electrical device with wet hands.

4. Products Management

4.3 If product Defect Occur

Situation	Confirmations and Solutions
No Power	<ul style="list-style-type: none">- Check the connection of power.- Check the Fuse-holder.
Increasing temperature in Bath	During the operation, temperature should be up to max. 70-80°C with no damages of equipment.
Off Heat-on indicator during Heat-on mode	Normal state (The time for sensor adjustment of the temperature to keep)
No Ultrasonic operation during Heat-on indicate	Caused the damages from Ultrasonic circuit and request A/S repair service.
※ If you have any problem other than the above, please call technical service of manufacturer or an official agent.	

5

5. Product Assurance

5.1 Quality Assurance

CERTIFICATE

of
**Quality Tested by Manufacturer
 and Free Warranty up to 2 years
 for Laboratory Instrument**

Limited Warranty

1. This instrument has gone through our testing process about all functions before shipment.
2. This instrument is warranted for 2 years from purchasing date.
3. A breakdown caused by customer's improper use or natural disaster is excepted from this warranty.
4. This certification is only for international customers.

Item	Analog Ultrasonic Cleaners
Model No.	WUC-A
Serial No.	
Date	
Customer	

DAIHAN Scientific Co., Ltd.

Daihan Bldg., 24-4, Sangwolgok-Dong
 Seongbuk-Gu, Seoul, KOREA, 136-120
 TEL: +82-2-967-5235, Ext. 2
 FAX: +82-2-963-5231
 Email: info@daihansci.co.kr
www.DAIHAN-Sci.com
www.ALLforLAB.com

Analog Ultrasonic Cleaners WUC-A

Certifications by ---

- ☑ CE Marked Products.
- ☑ ISO 9001 Certified.
- ☑ PL Insurance. (제조물 배상책임보험)

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